

UV/VIS Spectrophotometer

Basic Functions:

Model	Optical System	Wavelength Range	Bandwidth	Photo-metry	QTY	MULT	KINS	Scanning	Bio
BK-V1000	Single Beam	325-1020nm	4nm	•					
BK-UV1000	Single Beam	200-1020nm	4nm	•					
BK-V1200	Single Beam	325-1020nm	4nm	•	•				
BK-UV1200	Single Beam	200-1020nm	4nm	•	•				
BK-V1600	Single Beam	320-1100nm	4nm	•	•	•			
BK-UV1600	Single Beam	190-1100nm	4nm	•	•	•			
BK-V1800	Single Beam	320-1100nm	2nm	•	•	•	•		
BK-UV1800	Single Beam	190-1100nm	2nm	•	•	•	•		
BK-V1900	Single Beam	320-1100nm	2nm	•	•	•	•	•	•
BK-UV1900	Single Beam	190-1100nm	2nm	•	•	•	•	•	•
BK-S360	Single Beam	190-1100nm	1.8nm	•	•	•	•	•	•
BK-S380	Single Beam	190-1100nm	1nm	•	•	•	•	•	•
BK-S390	Single Beam	190-1100nm	0.5, 1, 2, 4, 5nm	•	•	•	•	•	•
BK-D560	Double Beam	190-1100nm	1.8nm	•	•	•	•	•	•
BK-D580	Double Beam	190-1100nm	1nm	•	•	•	•	•	•
BK-D590	Double Beam	190-1100nm	0.5, 1, 2, 4, 5nm	•	•	•	•	•	•

Features:

- * Large LCD screen (128*64 Dots).
- * Pre-aligned design ensures the user can change lamps conveniently.
- * With SiO₂ coating optical mirror, reducing the pollution from outside fully.
- * Environmental Deuterium lamp, preventing from Ozone inhalation.
- * High-class grating, wholly hermetic light path design, ensures the super low stray light.
- * Real-time monitoring the lifetime of Deuterium lamp and Tungsten lamp with advanced system.
- * Wavelength calibration, wavelength setting, change lamp source and dark current calibration automatically.
- * Optional PC Software to expand the applications to Quantitative, Multi-Wavelength and Kinetics.
- * Widely used for Organic chemistry, Inorganic chemistry, Life sciences, food, Medicine and health, Agriculture, Geology, Metallurgy and Environment etc. fields.



BK-UV/V1000/1200

Basic Functions:

* Photometry

Test Abs, Transmittance by fixed wavelength.

* Quantitative

- a. C Mode: Establish C=K*A curve, test sample concentration.
- b. F Mode: Establish A=K*(1/C+K0), test sample concentration.
- c. Display and save curve, test data (BK-V1200).

WL: 546.0 T: 100.0				WL: 546.0 C: 10C			
A	T	C	F	A	T	C	F

Photometry Mode

Con. Test C=K*A+B

Technical Parameters:

Model	BK-UV1000	BK-V1000	BK-UV1200	BK-V1200
Optical System	Single Beam, Grating 1200 lines/mm			
Wavelength Range	200~1100nm	325~1100nm	195~1100nm	320~1100nm
Spectral Bandwidth	4nm			
Wavelength Accuracy	±1.8nm		±1.0nm	
Wavelength Repeatability	±0.5nm			
Photometric Accuracy	±0.002A(0~0.5Abs), ±0.004A(0.5~1.0Abs), ±0.5%T(0~100%T)			
Photometric Repeatability	0.001Abs(0~0.5Abs), 0.002Abs(0.5~1.0Abs), ±0.2%T(0~100%T)			
Stray Light	≤0.2%T@360nm, 220nm(UV)			
Stability	±0.002A/h@500nm			
Noise	±0.001A			
Display	128*64 Dots LCD			
Photometric Mode	T, A, C, E			
Photometric Range	0~200%T, 0.301~3.0A			
Detector	Silicon Photodiode			
Light Source	Deuterium Lamp, Tungsten Lamp			
Input	Membrane Keypad			
Output	Parallel Print,USB1.0			
Power Supply	AC 85~265V, 50/60Hz			
Standard Accessories	1cm glass cuvette*4, 1cm quartz cuvette*2, 1cmquadruple cuvette holder*1	1cm glass cuvette*4, 1cmquadruple cuvette holder*1	1cm glass cuvette*4, 1cm quartz cuvette*2, 1cmquadruple cuvette holder*1	1cm glass cuvette*4, 1cmquadruple cuvette holder*1
	General application software, 1cm standard single slot cuvette holder, long optical path cuvette holder, cuvette, micro cuvette, test tube type absorber rack, solid sample holder, deuterium lamp(only for UV series), tungsten lamp, quadruple cuvette holder, printer, four-slot water bath type constant temperature absorption pool, X-Y Adjustable Micro Absorption Pool, single bath type constant temperature absorption pool			
Optional Accessories				
Package Size(W*D*H)	560*430*320mm			
Gross Weight	12kg	11.5kg	12kg	11.5kg